



ENVIKIT UREA



ENVIKIT 24Ure

From 50 mg/L

- Thank you for choosing products of Binh Lan ENVIBIOCHEM. Please read carefully before proceeding to ensure accurate results.
- Store in refrigerator.



1. Test directly on fish skin: press a test strip on wet fish, so that the strip became wet. Waite 2 minutes. The appearance of pink or pink-orange color is indication of urea.

☞ The pink or pink-orange color will disappear afterward.



2. Test of effluent from fish: Add 1.5 ml of effluent and 1 cm of test strip into Eppendorf tube. Waite 2 minutes. The appearance of pink or pink-orange color is indication of urea.

BINH LAN ENVIBIOCHEM COMPANY Limited

Web: <http://www.binhlan.com> **Email:** binhlanbiochem.gmail.com

Bank account: ACB Bank, Binh Lan Company, STK: 1246888

Office 1: Science-Technology Business Incubator, University of Science, HoChiMinh City
268 Ly Thuong Kiet Street, District 10, HoChiMinh City, Vietnam

Office 2: 595/208A Cach mang thang 8, Ward 15, District 10, HoChiMinh, Vietnam

Tel: 08-39700106, **Hotline:** 0985714580

For technical advice, price and ordering

Email to binhlanbiochem.gmail.com