



ENVIKIT FOR HARDNESS (GH)



ENVIKIT 03Hard 5 mg/l increment

- Thank you for choosing products of Binh Lan ENVIBIOCHEM. Please read instruction carefully before proceeding to ensure accurate results.
- Reagents contain strong base. In case of contact, immediately flush skin with plenty of water.



1. Fill the tube fully with sample water, then transfer water to the glass bottle.



2. Add 4 drops of Reagent 1. Swirl to mix.



3. Add 1 spoon of Reagent 2. Swirl to mix. Water turns purple.



4. Slowly add the Reagent 3 dropwise, swirling to mix after each drop while counting number of drops.



5. Continue dropping until water turns purple, then blue.



6. Stop when water turns blue.
GH (mg/l) = Number of drops × 10

- Wash all items by clean water, leave to dry in the air, and store in the box for the next test.
- If number of drops less < 4 then add another 1 tube of water to the glass bottle and continue steps 5 and 6.
GH (mg/l) = Number of drops × 10 : number of water tubes

BINH LAN ENVIBIOCHEM COMPANY Limited

Web: <http://www.binhlan.com> **Email:** binhlanbiochem.gmail.com

Bank account: ACB Bank, Binh Lan Company, 1246888

Office 1: Science-Technology Business Incubator of HoChiMinh City
268 Ly Thuong Kiet Street, District 10, HoChiMinh City, Vietnam

Office 2: 595/208A Cach mang thang 8, Ward 15, District 10, HoChiMinh, Vietnam

Tel: (84) -08-39700106, **Hotline:** (84)985714580

For technical support, quotation and order

Email to binhlanbiochem.gmail.com